



EasyGel: ECA Check® and ECA Check Plus®

ECA Check®:

- Supports growth and identification of three different groups of bacteria: E coli, coliforms and Aeromonas species all in one test.
- Isolation in 18-24 hrs
- Inhibits many Gram negative organisms and some Gram negative organisms

ECA Check Plus®

- Contains fluorescent compounds

Organism	Appearance- ECA Check	Appearance-ECA Check Plus
E coli	Dark blue to black with colored halo	Fluoresces bright blue
Coliforms	Light blue-gray to purplish pink	Does not fluoresce
Aeromonas	Pinkish red to light pink	Pinkish red to light pink
Salmonella	Light green	Does not fluoresce

ECA Check® and ECA Check Plus® as alternative to Coliscan® MF

- ECA Check® / ECA Check Plus® allow identification of 3 groups of bacteria simultaneously
- In order to separate Aeromonas from E coli using Coliscan® MF you must use specific inhibitors-antibiotic or high incubation temperature

John R Beck PhD
Weber Scientific
May 29, 2007